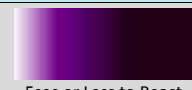
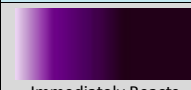

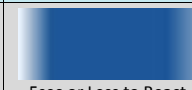
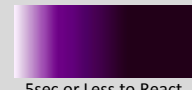
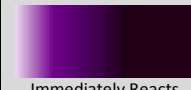
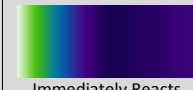


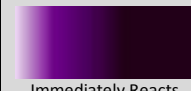
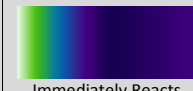
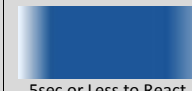
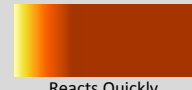
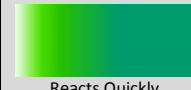
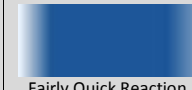
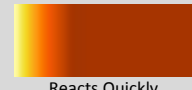
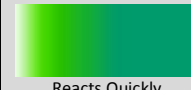




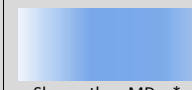

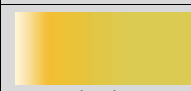
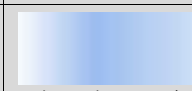
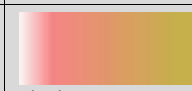






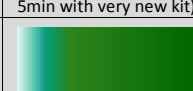
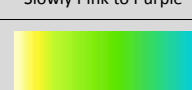
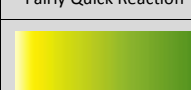
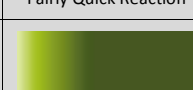

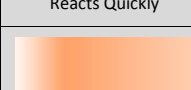

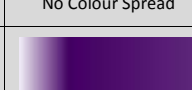

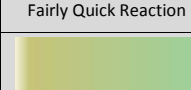
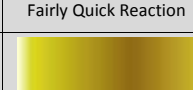
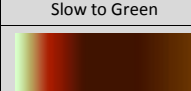
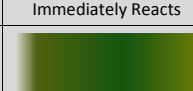

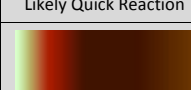
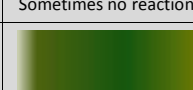
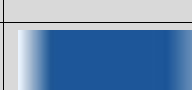
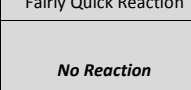


MARQUIS	MANDELIN	MECKE	SIMON'S	ROBATEST	SUSPECTED DRUG	NOTES
 5sec or Less to React	 Immediately Reacts	 Immediately Reacts	 5sec or Less to React	<i>No Reaction</i>	<b>MDMA or MDEA</b>	>10sec to react with Marquis or Mandelin indicates very low potency.
 5sec or Less to React	 Immediately Reacts	 Immediately Reacts	<i>No Reaction</i>	 Reaction Takes 1-5mins	<b>MDA</b>	
 Immediately Reacts	 Immediately Reacts	 Immediately Reacts	 5sec or Less to React	<i>No Reaction</i>	<b>MDMA or MDEA &amp; Methamphetamine</b>	May need to test a couple of times with Marquis to see yellow/orange reaction.
 Reacts Quickly	 Reacts Quickly	<i>No Reaction</i>	 Fairly Quick Reaction	<i>No Reaction</i>	<b>Methamphetamine</b>	
 Reacts Quickly	 Reacts Quickly	<i>No Reaction</i>	<i>No Reaction</i>	 Reaction Takes 1-5mins	<b>Amphetamine</b>	
 Immediately Reacts	 Immediately Reacts	 Slowly Changes Colour	 Slower than MDxx*	 Small Brown Ring	<b>Methylone (bk-MDMA)</b>	
<i>No Reaction</i> (Some Impure Batches React Yellow)	 Immediately Reacts	<i>No Reaction</i> (may smoke)	 Slower than MDxx*	 Slowly to a Brown Ring	<b>Mephedrone (4-MMC)</b>	
 Reacts Quickly	 Immediately Reacts	 Immediately Reacts	<i>Unknown</i>	<i>Unknown</i>	<b>LSD</b>	Best results seen testing LSD liquid (drops). If testing a pill use a large scraping. Won't test LSD tabs (blotters).
<i>No Reaction</i>	 No Reaction with 1yr Old Kit! USE NEW KIT!	<i>No Reaction</i> (Pale light yellow after 5min with very new kit)	<i>No Reaction</i>	<i>No Reaction</i>	<b>Ketamine</b>	Sometimes PMA shows no reaction to Mecke and could be mistaken for Ketamine!
 Slowly Pink to Purple	 Fairly Quick Reaction	 Fairly Quick Reaction	<i>No Reaction</i>	<i>No Reaction</i>	<b>Heroin</b>	If not much powder is used, Mandelin reaction can start as thin red brown streaks slowly changing to brown.
 Blue after a few mins	 Reacts Quickly	 Reacts Quickly	<i>No Reaction</i>	 No Colour Spread	<b>2C-B</b>	Check Acinad Romanoff channel on YouTube for reactions to other 2C-x or 2C-T-x chemicals
<i>No Reaction</i>	 Fairly Quick Reaction	 Fairly Quick Reaction	<i>No Reaction</i>	 No Colour Spread	<b>2C-T-7</b>	2C-T-7 -dangerous if consumed with MDxx* or a number of other drugs
 >10sec to React	 Slow to Green	 Immediately Reacts	<i>No Reaction</i>	<i>No Reaction</i>	<b>DXM</b>	DXM -dangerous if consumed with MDxx* or a number of other drugs
<i>No Reaction</i>	 Likely Quick Reaction	 Sometimes no reaction	<i>No Reaction</i>	 Expected slow Reaction	<b>PMA</b>	PMA - high risk of death if consumed
<i>No Reaction</i>	 Fairly Quick Reaction	 Reacts Quickly	 Fairly Quick Reaction	<i>No Reaction</i>	<b>PMMA</b>	PMMA - high risk of death if consumed
<i>No Reaction</i>	<i>No Reaction</i>	<i>No Reaction</i>	<i>No Reaction</i> (Or May Have a Faint Blue Reaction)	<i>No Reaction</i>	<b>Piperazines (such as BZP / TFMP / mCPP)</b>	Piperazines - high risk of bad side effects if consumed and some risk of death
<i>No Reaction</i>		<i>No Reaction</i>	<i>Unknown</i>	<i>Unknown</i>	<b>Strychnine</b>	Very toxic poison - high risk of death if consumed

EVEN USING THIS CHART SOME REACTIONS MAY LOOK THE SAME FOR SOME DIFFERENT MIXTURES OF DRUGS AND OR IMPURITIES!

(Version: 18-Dec-2019)

Example: MDxx + PMA or PMMA – may appear same reaction as containing only MDxx. Larger risk of death than consuming only PMA or PMMA!

Example: MDxx + Piperazines – will appear same reaction as containing only MDxx. High risk of bad side effects if consumed and some risk of death.

To watch videos of many different, lab verified drug samples reacting to the different reagents please check out: [Acinad Romanoff](#) channel on YouTube.

Please check [pillreports.net](#) and [ecstasydata.org](#) to see if there are any reports on your pill.